


Assigning the detector to channel


FV3000

- 1 Press the **Dye & Detector Select** button on [PMT Setting] Tool Window. The [Dye & Detector Select] dialog box appears.

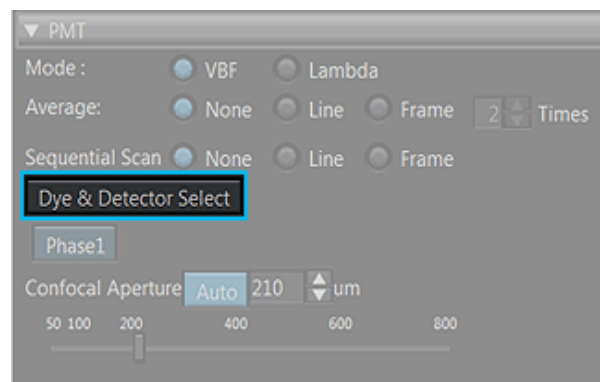


- 2 Select the [Dye] or the [Channel] tab in the [Dye & Detector Select] dialog box.


| | |
|---------|---|
| Dye | Specify the fluorescent reagent (Dye) and assign the observation channel. |
| Channel | Specify the detector and assign the observation channel. |

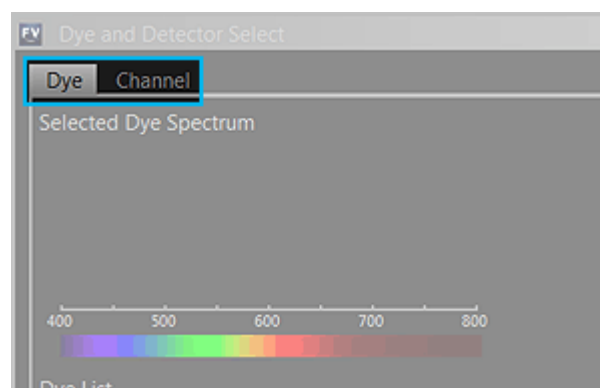


- 3 If you want to observe in multiple phases using the virtual channel scan, press the **Add Phase** button.
(The virtual channel scan is the option of FV3000.)

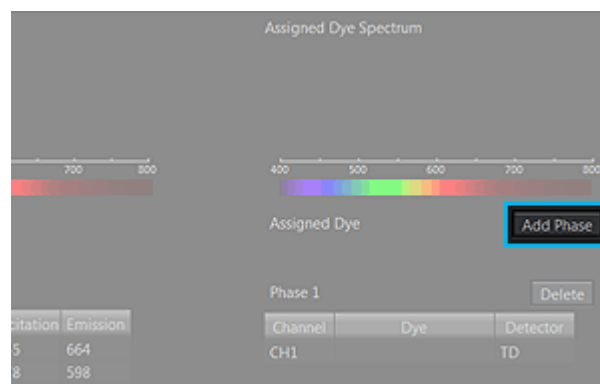


- 4 If you selected [Channel] in
 - 2, select the detector to be assigned to the observation channel from [Detector List].

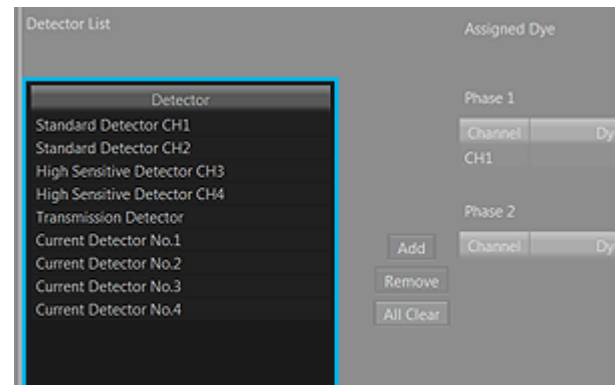

If you selected [Dye] in **2**, select "Dye" to be observed from [Dye List].

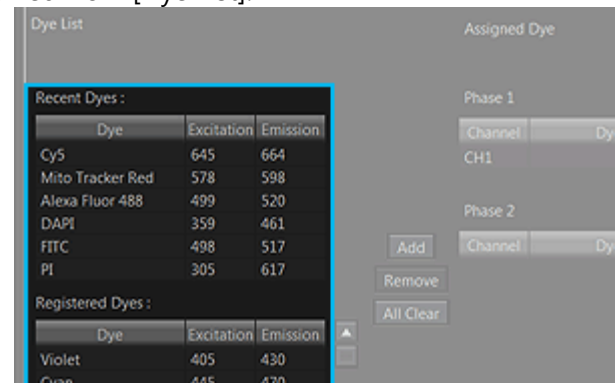

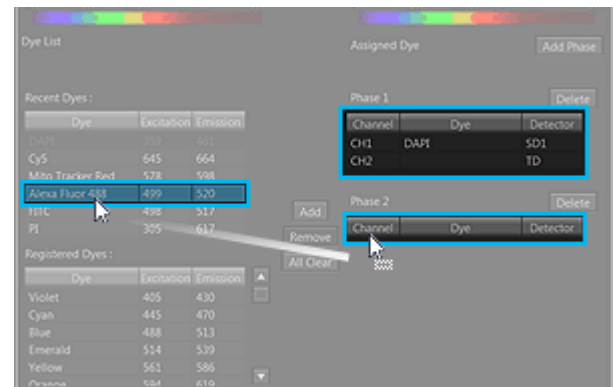

The observation channel number and the abbreviation (physical channel name) of the assigned detector are displayed in the observation channel list.

6

After setting all channels, press the **OK** button.

