



## Recording 3D image movement

The movie can be created by moving the 3D image by dragging the mouse.


- 1 Select [Volume Setting] in the [Tool Window] menu. [Volume Setting] Tool Window is displayed.

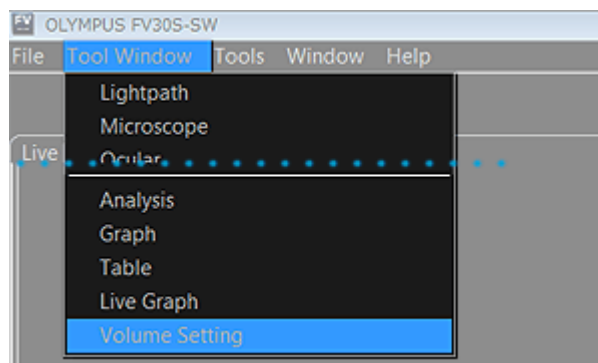


- 2 Select [\[View\] tab in \[Volume Setting\] Tool Window](#).  
(2 and 3 are the procedures necessary for moving the 3D image by dragging the mouse.)

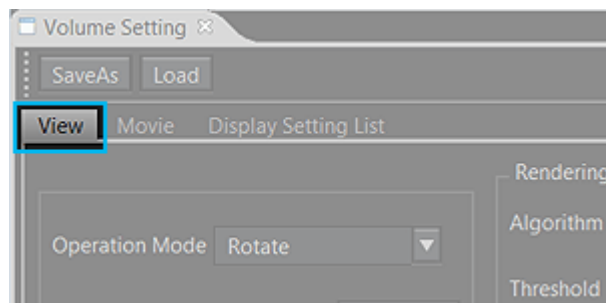

- 3 Select "Rotate" in [Operation Mode].

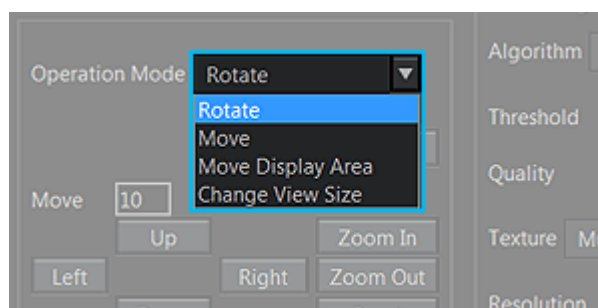


- 4 Select [\[Movie\] tab in \[Volume Setting\] Tool Window](#).

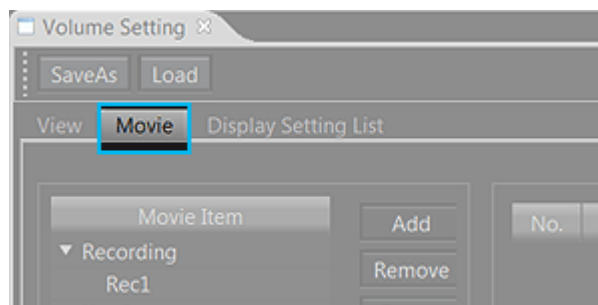


5


Select "Recording" in [Movie Item] and press the **Add** button. "Rec1" is shown below "Recording".

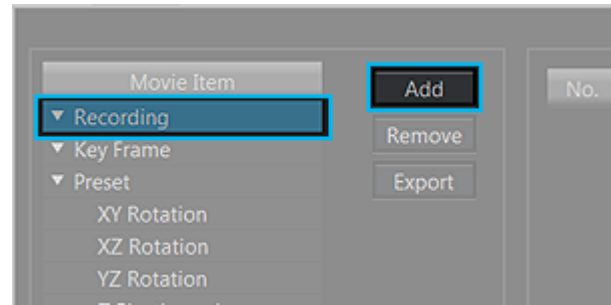


**6** Select "Rec1" in [Movie Item].

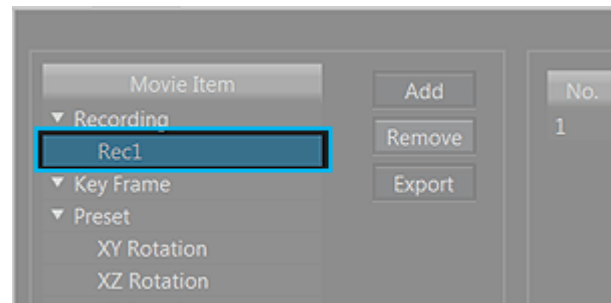


**7** Double-click on the value of [FrameRate] to display the ▼ button. Press the ▼ button to select the frame rate.

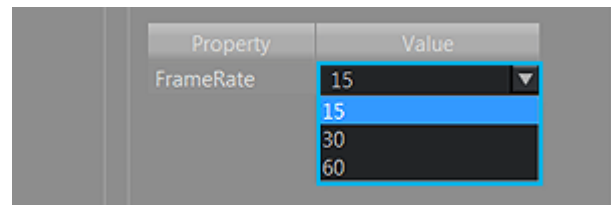

**8** Press the **Rec** button to start recording, and move the 3D image in [\[Image\] window](#) by dragging.

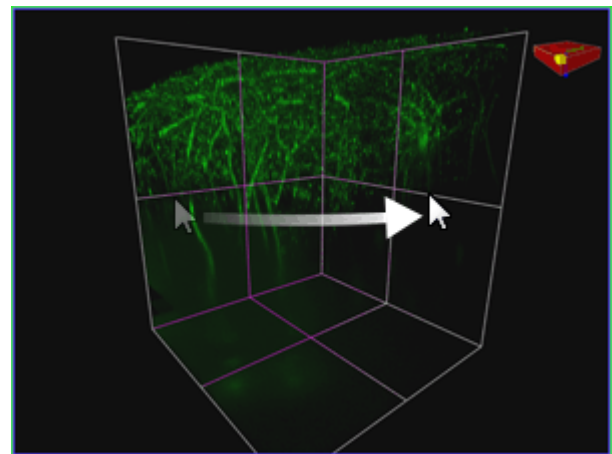

**9** Press the **Stop** button to stop recording.

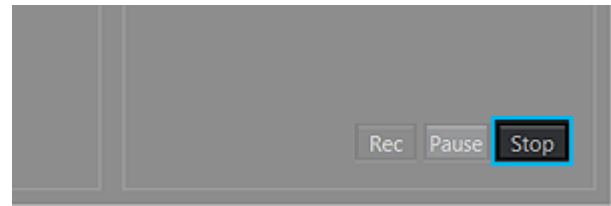**10**

Press the **Play** button to play back the movie.

