



Exporting movie


The movie can be exported (saved) in WMV type.

- 1 Select "Rec xx" or "Sequence xx" in [Movie Item] and select the movie you want to export.
 Rec xx : This is a movie recorded with moving the 3D image by dragging the mouse.
 Sequence xx : This is a movie created by interpolating between Key Frames after Key Frames are registered with moving the 3D image by dragging the mouse.

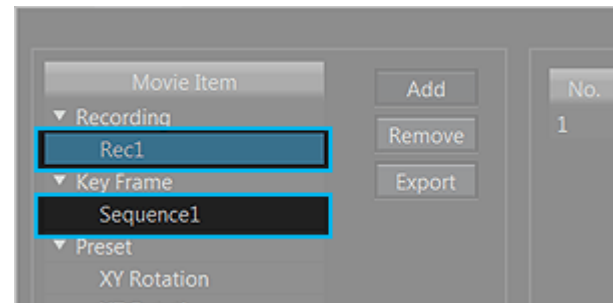


- 2 Press the **Export** button. The [Export] dialog box appears.

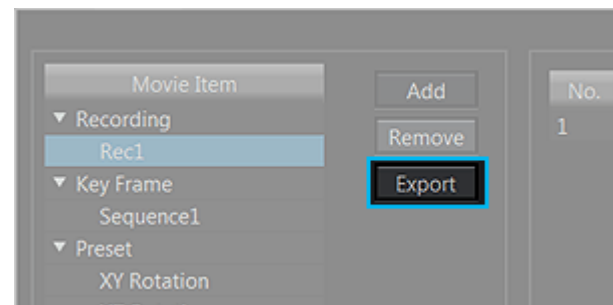


- 3 The save destination of the movie to be exported is displayed in **a**. If you want to change it, press the  button to display the dialog box and select the folder of the save destination.

 In order to organize movie files with ease, it is recommended to specify the save destination.

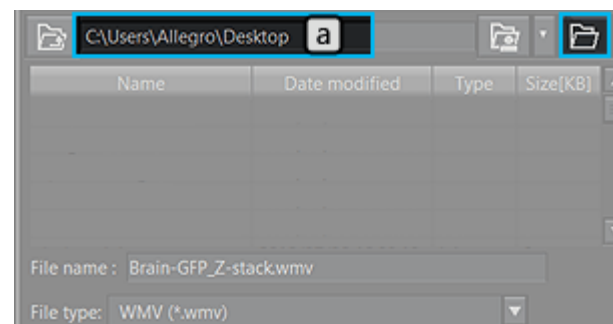


- 4


The file name of the movie to be exported is displayed in [File name]. If you want to change it, enter the file name directly.

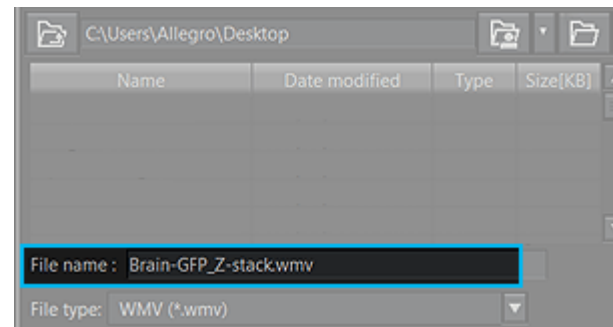

- 5 The default value is displayed in [Frame rate]. If you want to change it, enter the value directly.

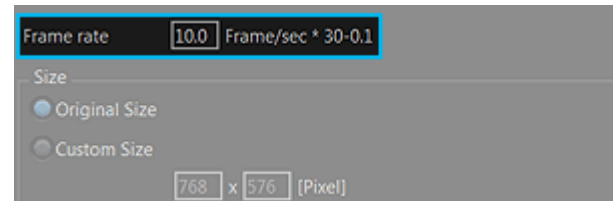

- 6 If you export the image by changing the original image size, select "Custom Size" in [Size] and enter the value directly.

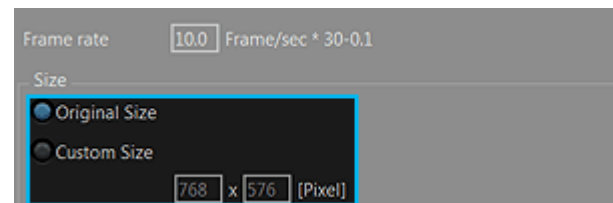

- 7 Press the **Save** button. The movie is exported.

